6 | GLASS FILTRATION APPARATUS AND ACCESSORIES

Due to their high chemical and thermal shock resistance DURAN® filters and the corresponding filter plates are ideal for separations, e. g. with strong acids or alkalis. Thus they offer advantages in comparison with other materials such as plastic or paper. DURAN® filter products have a maximum operating temperature of +450 °C and are therefore far superior to other materials.

The corresponding filtration vessels are specially optimised to the matching filtration apparatus and are vacuum-tight due to their special geometry and high wall thickness. This characteristic has been approved by the German TÜV and marked with the "GS"-indication.

The glass filters are classified as being in porosity classes 0 to 5 according to their nominal maximum pore size. The following table shows the corresponding porosity range. The specified pore sizes always relate to the largest pore in the plate. This specification also characterises the minimum nominal size of particles which may be retained by the filtration.

Usage tips:

- The maximum permissible operating temperature is +450 °C.

- Uniform heating is recommended to avoid thermal stresses and resultant breakages.
- Heat glass filtration apparatus with disk diameters of more than 20 mm in initially cold ovens or sterilisers only.
- The heating or cooling rate should not exceed 8 °C/min.
- When filtering hot substances observe the thermal shock resistance and, if necessary, preheat the filtration apparatus in a drying cabinet.
- Wet filtration apparatus should be heated slowly up to +80 °C and dried for one hour before increasing the temperature further.

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Particular suited to use with hot or aggressive media thanks to the thermal shock and chemical resistance of DURAN®. Also available in **soda-lime glass²**.

Typical applications: decanting and filtration of substances.

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From DURAN® glass with its good thermal shock and chemical resistance. Also available in **soda-lime glass¹**.

Typical applications: decanting of powdered substances and granulated material.

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
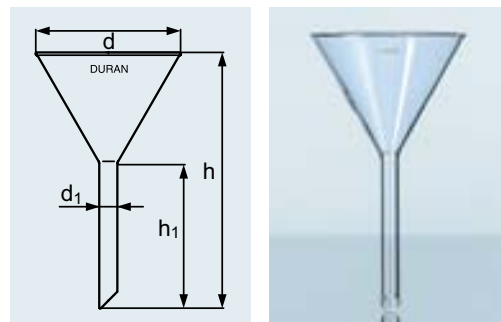
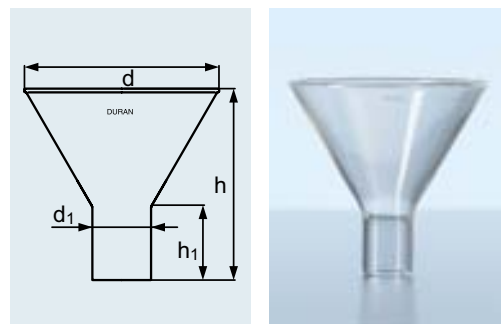
DURAN® POWDER FUNNEL with short, wide stem

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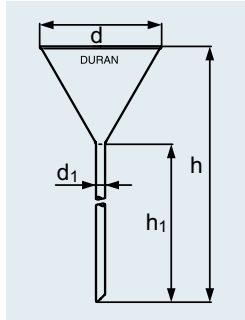From DURAN® glass with its good thermal shock and chemical resistance. Also available in **soda-lime glass³**.

Typical applications: filtering and decanting of liquids of different densities.

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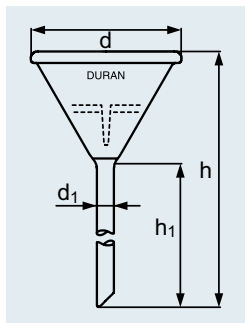
¹ Non-DIN ISO size.

² Details apply only to DURAN® version.

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DURAN® ANALYTICAL FUNNEL
 for quick filtration

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From DURAN® glass with its good thermal shock and chemical resistance.

Typical application: for rapid liquid filtration.

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From DURAN® glass with its good thermal shock and chemical resistance. The ribbed form is ideal for filtering with round-paper.

Typical application: filtering of liquids.

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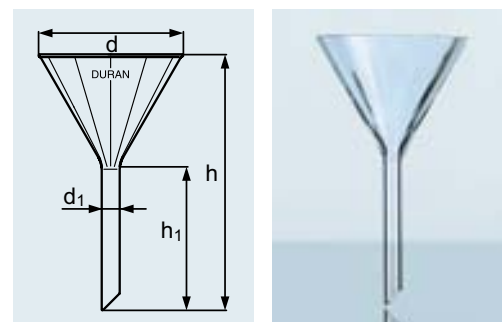
Made from DURAN® glass with its good thermal shock and chemical resistance. Filter funnels mate to the filtering flask via a conical rubber seal (GUKO).

Typical application: filtering of liquids.

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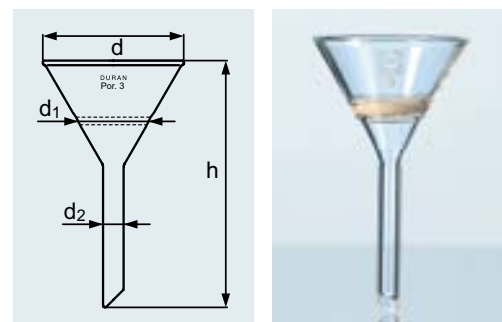
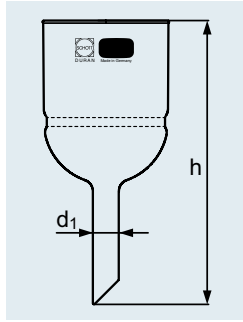
From DURAN® glass with its good thermal shock and chemical resistance. Filter funnels mate to the filtering flask via a conical rubber seal (GUKO).

Typical applications: qualitative inorganic analysis and preparative chemistry.

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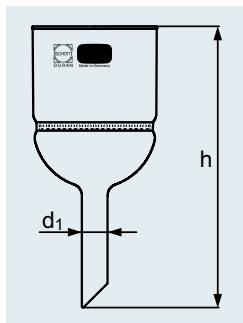
From DURAN® glass with its good thermal shock and chemical resistance. The Buechner funnel features a glass support for membrane and paper filters.

Best.-Nr. Cat. No.	Inhalt Capacity ml	Platten Disc d mm	d _i mm	h mm	passende Papier- Rundfilter matching filter paper d mm	AD OD mm	Ver- packungs- einheit/Stück Pack/Quantity
21 341 22	70	48	10	132	45	57	1
21 341 28	125	60	10	140	55	72	1
21 341 34	220	73	18	190	70	90	1
21 341 44	500	95	22	240	90	106	1
21 341 54	1 000	120	22	270	110	136	1

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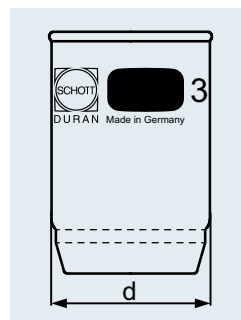
¹ Substitute required porosity (see page 239) for X in the Cat No.


Best.-Nr. Cat. No.	d mm	passender Filtriervorstoß Best.-Nr. suitable filter adapter Cat. No.	Verpackungseinheit/Stück Pack/Quantity
29 201 14	26	24 316 16	10
29 201 21	33	24 316 22	10
29 201 26	41	24 316 26	10
29 201 31	49	24 316 32	10

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¹ For chemical and thermal resistance, see page 222.

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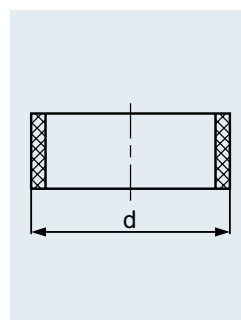
RUBBER SLEEVE from EPDM¹ for filter crucibles

Best.-Nr. Cat. No.	d mm	d _i mm	h mm	passende Gummi- manschette Best.-Nr. suitable rubber sleeve Cat. No.	Verpackungseinheit/ Stück Pack/Quantity
24 316 16	27	10	108	29 201 14	10
24 316 22	34	10	110	29 201 21	10
24 316 26	41	10	125	29 201 26	10
24 316 32	50	10	132	29 201 31	10

DURAN® FILTRIERVORSTOSS

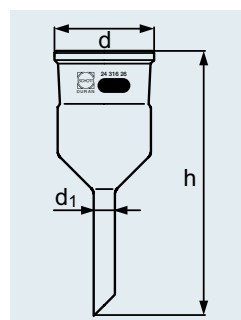
DURAN® FILTER CRUCIBLE / FUNNEL ADAPTER

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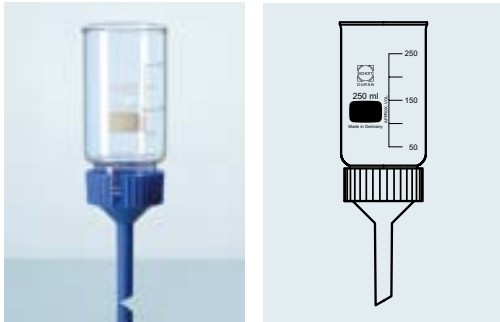**DURAN® FILTER FUNNEL HEAD
 WITH PP¹ FUNNEL AND TWO
 FKM¹ SEALS**

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Interchangeable filter disks. Available in three filter diameters, each of which is available in four different porosities. Important: insert filter disc between two FKM seals. After filtration, disk can be removed to allow simple and safe removal of the filtrand. Long filter disk service life, as disks are not damaged when the filtrand is scraped off. Easy cleaning of both sides is possible. Cost-effective as components and disks can be ordered individually as required.

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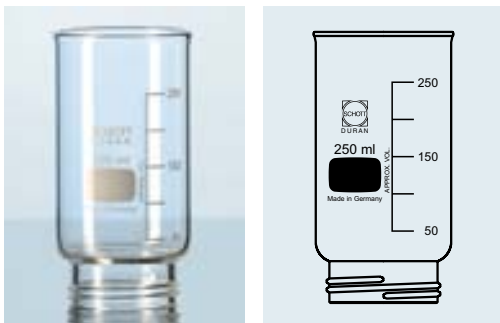
From DURAN® glass with its good thermal shock and chemical resistance. Fused glass rim.

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
DURAN® FILTER DISK WITH GLASS RIM
patented

Best.-Nr. Cat. No.	Porosität Porosity	für Platte for disc d mm	Verpackungseinheit/ Stück Pack/Quantity
25 202 4X ¹	1-4	24	1
25 205 0X ¹	1-4	50	1
25 209 0X ¹	1-4	90	1

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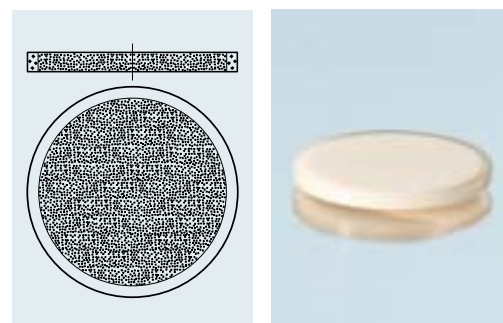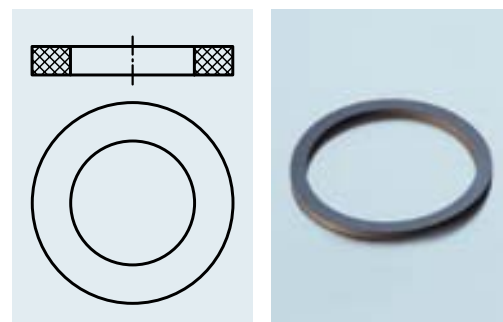
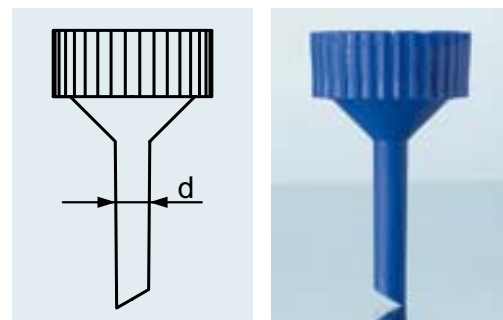
FUNNEL FOR FILTER FUNNEL HEAD from PP¹

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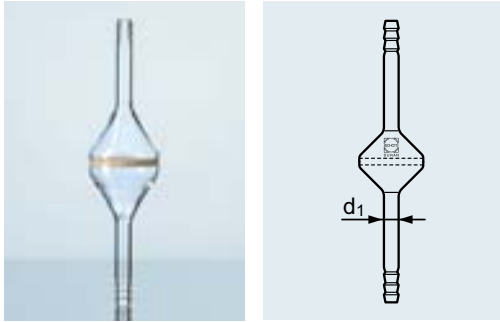
Typical application: in-line filtration of gas lines to remove solid impurities (e.g. dust).

Best.-Nr. Cat. No.	Porosität Porosity	Platte Disc d mm	d ₁ mm	Verpackungseinheit/Stück Pack/Quantity
25 855 0X ¹	I-4	30	10	I
25 855 1X ¹	I-4	60	15	I
25 855 2X ¹	I-4	90	16	I

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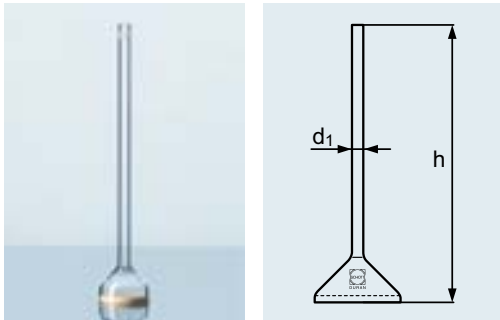
DURAN® IMMERSION FILTER
for reverse filtration

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Typical application: extraction of clear filtrate (not the filtrand).

Best.-Nr. Cat. No.	Porosität Porosity	Platte Disc d mm	d ₁ mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 855 6X ¹	I-4	35	10	210	I

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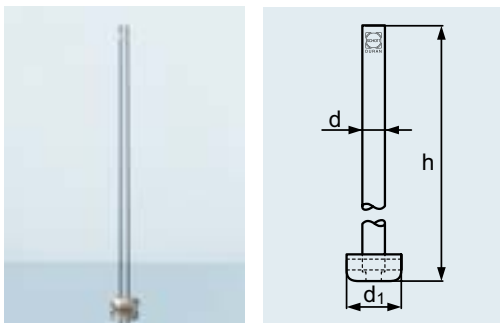
Typical application: reactions between gases and liquids.

Best.-Nr. Cat. No.	Porosität Porosity	d mm	d ₁ mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 856 0X ¹	0-2	6	22	250	5
25 856 1X ¹	0-2	9	25	250	5
25 856 2X ¹	I	11	34	250	5

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With screw-connection system. The insertion height of the head is adjustable. Individual parts can also be ordered separately (see page 140).

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
Typical applications: cleaning (“washing”) of gases with solvents.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	Olive Hose connection d mm	Verpackungseinheit/ Stück Pack/Quantity
24 713 00	500	45	9	1

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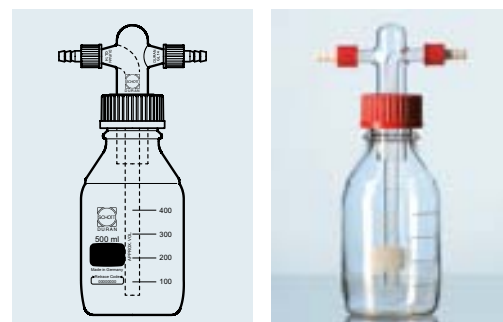DURAN® GAS WASHING BOTTLE
head Drechsel-type, without filter disk

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With screw-connection system. The insertion height of the head is adjustable. The distribution of the gas in the liquid is carried out through the filter disk and thus the absorption efficiency significantly increased. Gas washing bottles also work reliably at high flow rates. Individual parts can also be ordered separately (see page 140).

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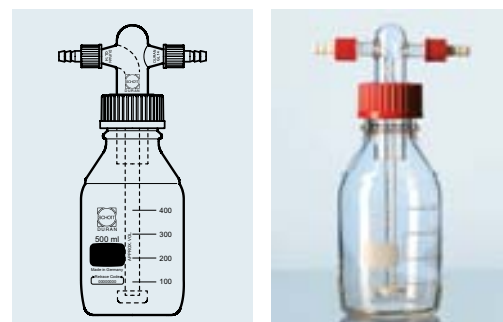
DURAN® GAS WASHING BOTTLE
head Drechsel-type, with filter disk

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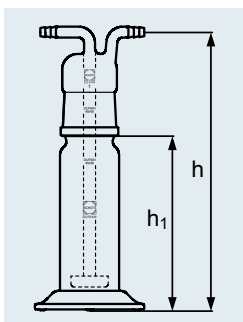
DURAN® GAS WASHING BOTTLE
head with filter disk, with standard ground joint

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
Typical application: cleaning („washing“) of gases with solvents.

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DURAN® GAS WASHING BOTTLE
 with fused-in filter disk,
 with standard ground joint and cap

Best.-Nr. Cat. No.	Porosität Porosity	Inhalt Capacity	NS Joint	d	h	h ₁	Olive Hose connection d mm	Verpackungs- einheit/Stück Pack/Quantity
25 703 01	I	350 ml	29/32	60 mm	250 mm	180 mm	10	I

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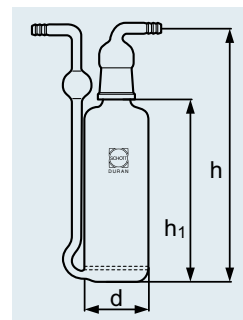
DURAN® MICRO IMMERSION FILTER
 for reverse filtration

Best.-Nr. Cat. No.	Porosität Porosity	d ₁ mm	d mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 857 1X ¹	I-4	10	6	100	10

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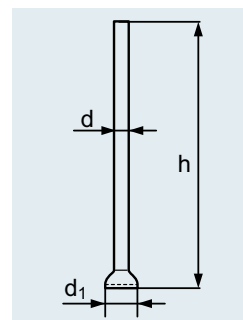
¹ Substitute required porosity (see page 239) for X in the Cat. No.


DURAN® MIKRO-FILTERKERZE
 ohne Rohr

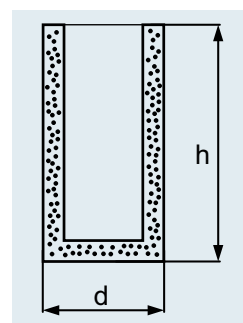DURAN® MICRO FILTER CANDLE
 without tube

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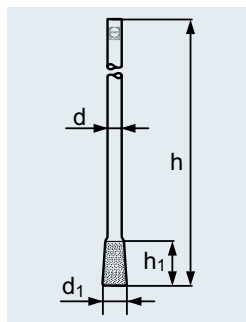DURAN® MICRO FILTER CANDLE
with narrow tube

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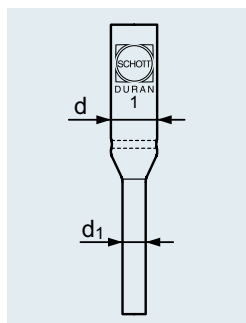
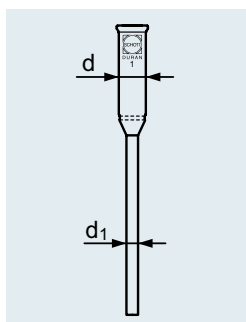
DURAN® MICRO FILTER FUNNEL
Pregl type

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
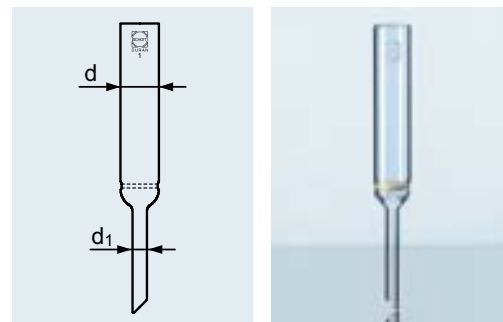
Heavy walled for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". For suitable rubber gaskets, see page 147.

Typical applications: separations by vacuum filtration.

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DURAN® FILTERING FLASK WITH SIDE-ARM SOCKET
Erlenmeyer shape

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	d ₁ mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
21 183 36	250	85	34	155	10
21 183 44	500	105	34	185	10
21 183 54	1 000	135	45	230	10
21 183 63	2 000	166	60	255	1

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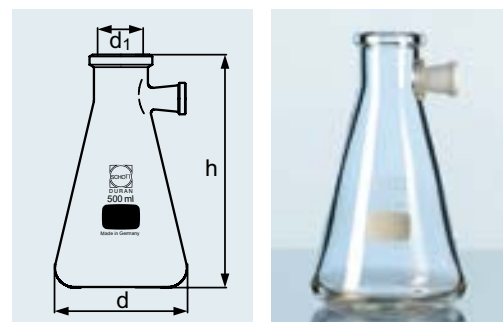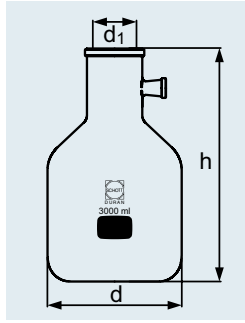
DURAN® FILTERING FLASK WITH SIDE-ARM SOCKET
bottle shape

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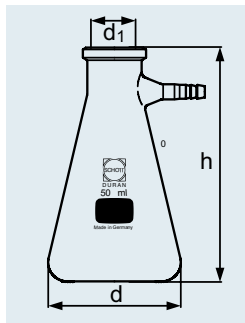
Note: These filtering flasks have a ground socket 17.5/26 for vacuum tube of 15 to 18 mm OD (e.g. 6 x 5 mm or 8 x 5 mm, DIN 12 865).

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Provision of filtration flasks with a socket has not only made work in preparation and analytical laboratories easier and simpler, but has also reduced the risk of accidents.

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DURAN® FILTERING FLASK WITH GLASS HOSE CONNECTION
Erlenmeyer shape

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Due to the heavy wall thickness the apparatus is vacuum-tight. For suitable rubber gaskets, see page 147.

Typical application: separations by vacuum filtration.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	d ₁ mm	h mm	Olive Hose connection d mm	Verpackungseinheit/ Stück Pack/Quantity
21 201 24 ¹	100	64	24	105	11	10
21 201 36 ¹	250	85	34	155	11	10
21 201 44 ¹	500	105	34	185	11	10
21 201 54 ¹	1 000	135	45	230	11	10
21 201 63 ¹	2 000	166	60	255	11	1

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Heavy walled for vacuum use. For suitable rubber gaskets, see page 147.

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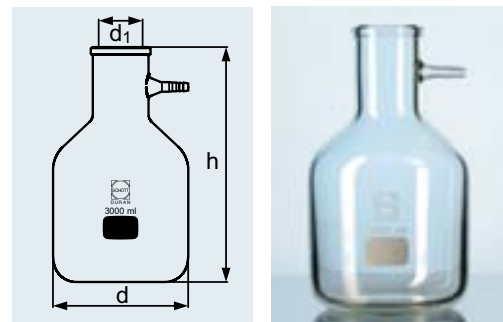
¹ Does not conform to the "equipment and product safety regulations".

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Flaschenform

DURAN® FILTERING FLASK WITH GLASS HOSE CONNECTION bottle shape

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Heavy walled for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". The plastic hose connection is replaceable. For suitable rubber seals, see page 147.

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
Typical application: Separations by vacuum filtration.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	d ₁ mm	h mm	Olive Hose connection d mm	Verpackungs- einheit/Stück Pack/Quantity
21 204 24 5	100	64	24	105	9	10
21 204 36 5	250	85	34	155	9	10
21 204 44 5	500	105	34	185	9	10
21 204 54 5	1 000	135	45	230	9	10
21 204 63 5	2 000	166	60	255	9	1

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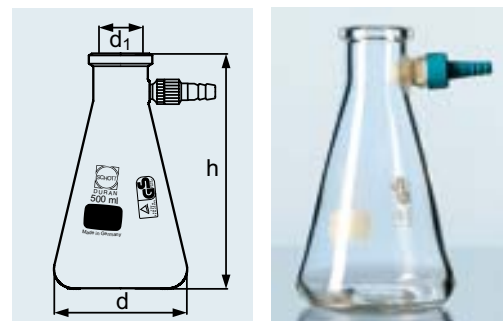
Flaschenform

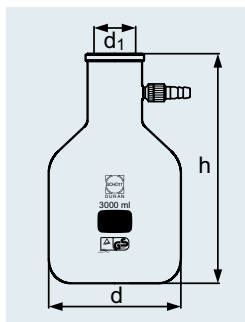**DURAN® FILTERING FLASK
 WITH KECK™ ASSEMBLY SET**
bottle shape

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Due to the heavy wall thickness the apparatus is for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". The plastic hose connections can be replaced. For suitable rubber gaskets, see page 147.


Typical application: separations by vacuum filtration.

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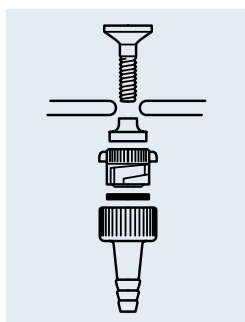*With removable plastic hose connection (PBT¹), short and long screw (PP¹), seals (VMQ¹, EPDM¹).
 Suitable for filtering flasks 100 - 20 000 ml.*

Best.-Nr. Cat. No.	Olive Hose connection d mm	Verpackungseinheit/Stück Pack/Quantity
29 258 54	9	10

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
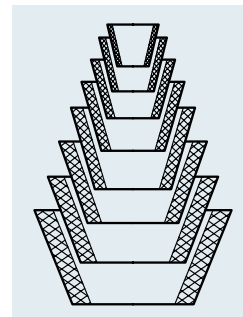
GUKO FROM EPDM¹
conical rubber gaskets for filtering flasks

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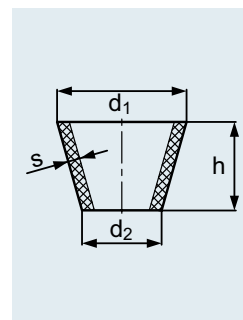
Virtually universal applications, as the medium only comes into contact with glass and PTFE. The scaled funnel simplifies dosing and analysis. With PTFE plate holder. Filter paper, membrane filters (47 mm) or glass filters can be used for filtration. Plates and PTFE adapters are replaceable. Easy and fast cleaning. All components are available as spare parts.

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Application examples: Coarse and fine filtration, filtration of HPLC media, residue analysis

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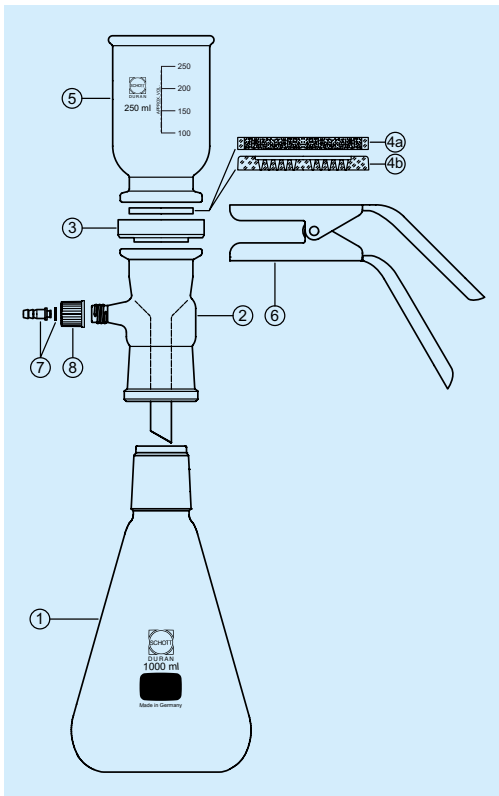
¹ Comprises the following order numbers:
 21 204 445, 29 202 27, 24 316 26, 29 201 26,
 25 851 23 and 25 851 24.

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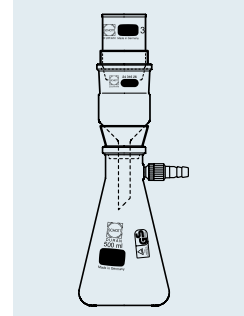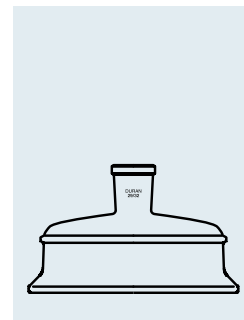
DURAN® LID FOR FILTER
APPARATUS, WITT TYPE
with tubulature, standard ground joint 29/32

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*DURAN® FILTER APPARATUS,
 WITT TYPE*

*complete with interchangeable lid and KECK™
 assembly set¹, standard ground joint 29/32*

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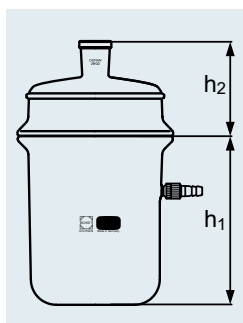
*DURAN® FILTER APPARATUS,
 WITT TYPE*

*complete with interchangeable lid and KECK™
 assembly set¹, standard ground joint 45/40*

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
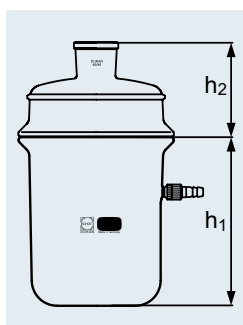
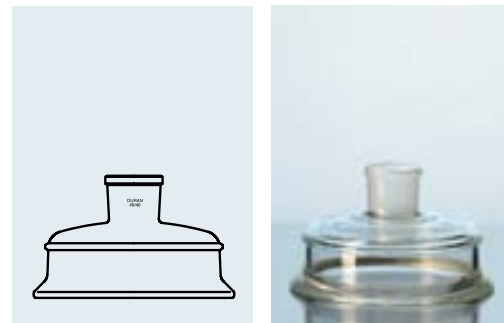
*DURAN® LID FOR FILTER APPARATUS,
WITT TYPE*
*wide rough-ground tubulature, standard
ground joint 45/40*

Best.-Nr. Cat. No.	Wasserverbrauch mind. Water flow rate min. l/h	Wasserdruck mind. Water pressure min. bar	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 362 99	300	1,2	275	1

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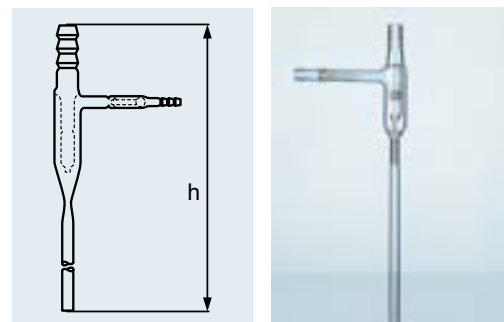*DURAN® WATER JET PUMP
with non-return valve*

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*Throughput: 400 l/h at 3.5 bar water
pressure und 12 °C water temperature.*

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*WATER JET PUMP
from plastic (PP¹), with non-return valve,
hose connection and adapters for 1/2" and 3/4"*

Best.-Nr. Cat. No.	Olive passend für Schlauch ID Connection suitable for hose ID mm	Wasserverbrauch mind. Water flow rate min. l/h	Wasserverbrauch mind. Water flow rate min. bar	h mm	Verpackungs- einheit/Stück Pack/Quantity
29 250 01	9-12	170	1	235	1

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